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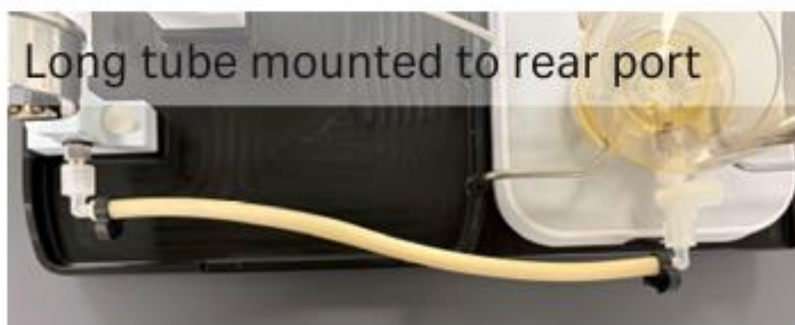
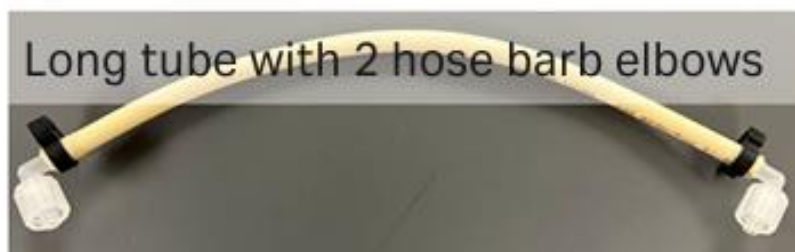
Cytiva
100 Results Way
Marlborough, MA 01752
United States of America

Notification date: May 14, 2026

Change notice: Minimate™ EVO TFF System 3-way valve change

Dear valued customer,

A discontinuation of the original 3-way valves used in the Minimate TFF System has prompted a replacement. Effective immediately, Minimate systems will include 3-way valves of the same material as the original, but with a mirror image orientation of port 2 as laid out in our document: *System setup options with updated 3-way valves*. In this document we explain multiple ways to account for the new valves during system setup. Recommended tubing lengths and additional connectors are described in detail with photo examples.

We will be adding an additional male luer lock 1/8" hose barb elbow to our Minimate fittings kit and to accompany new system purchases. We'll be working to update our product literature on the web to include the new instructions but, in the immediate term, please use the *System setup options* document referenced in this letter to ensure correct setup.

We apologize for any confusion this has caused. Please contact your local Cytiva sales team for any questions related to this change. Thank you for your continued support and business with Cytiva.

Sincerely,

A handwritten signature in black ink, appearing to read 'Joshua Smith', written in a cursive style.

Joshua Smith
Global Product Manager – Cellular
Cytiva
joshua.smith@cytiva.com

© Copyright 2026 Cytiva. Cytiva and the Drop logo are trademarks of Life Sciences IP Holdings Corporation or an affiliate doing business as Cytiva. Minimate is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva. **This Notice is proprietary to Cytiva and is furnished in confidence for the private use of the intended recipient, for the sole purpose of evaluating Cytiva technology. It may not be copied nor disclosed, in whole or in part, without the prior express approval of Cytiva.**

cytiva.com

Confidential - Company Proprietary

Minimate™ EVO TFF system

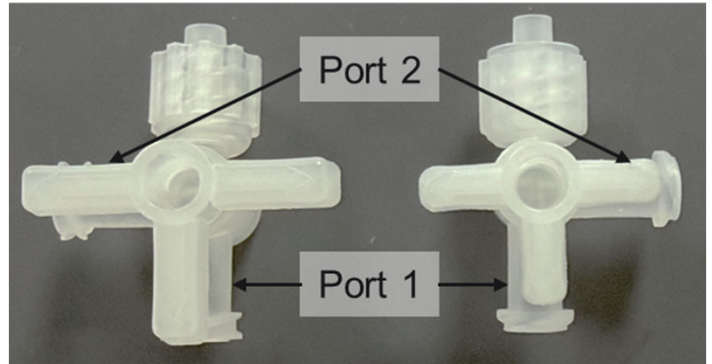
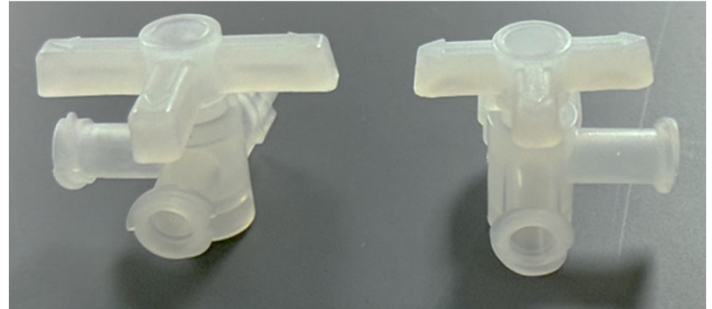
System setup options with updated 3-way valves

Introduction

Discontinuation of the 3-way valve (Old) used with the Minimate EVO TFF system led to its replacement with a similar type valve (New) that is its mirror image with respect to the port 2 outlet.

The changed port orientation has consequences for instrument setup. This document describes ways to workaround the current geometry.

One solution for the valve at the front reservoir port and two solutions for the valve at the rear reservoir port.

Old

New


Use of 3-way valves on Minimate EVO TFF system

The 3-way valves are used at the front and rear ports of the reservoir.


Rear port – The long tube is connected to port 2

Front port – The short tube is connected to port 1

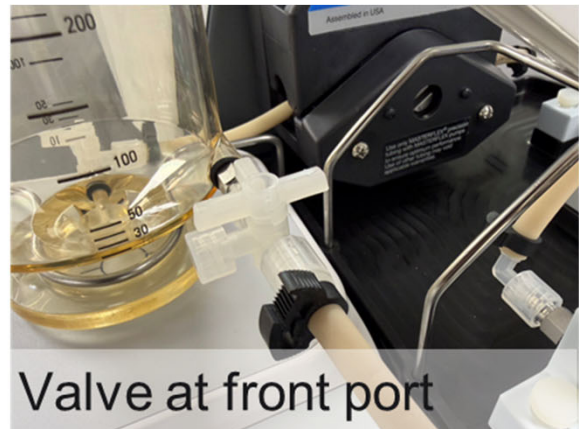Connections here are shown with the discontinued 3-way valves

Drip tray assembly naming conventions

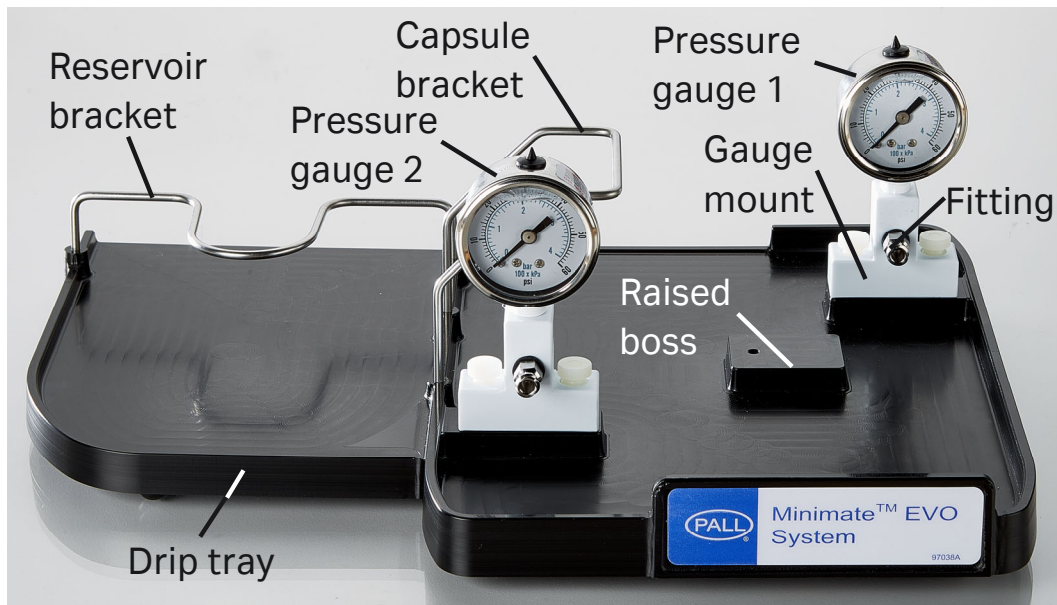The drip tray has three raised bosses, two of which are occupied with pressure gauges that are attached via gauge mounts.

Fittings on the mount are used to connect tubing (not shown).

The third boss is used to attach the capsule stand (not shown)

The reservoir and capsule brackets complete the drip tray assembly.

Tubing lengths for Options 1 and 2

Tubing	Length	
	Option 1	Option 2
Long Tube	8½" (23.5 cm)	9¼" (21.6 cm)
Mid-length tube	6" (15.2 cm)	6" (15.2 cm)
Short tube	3" (7.6 cm)	3" (7.6 cm)

Recommendation for 3-way valve at reservoir front port

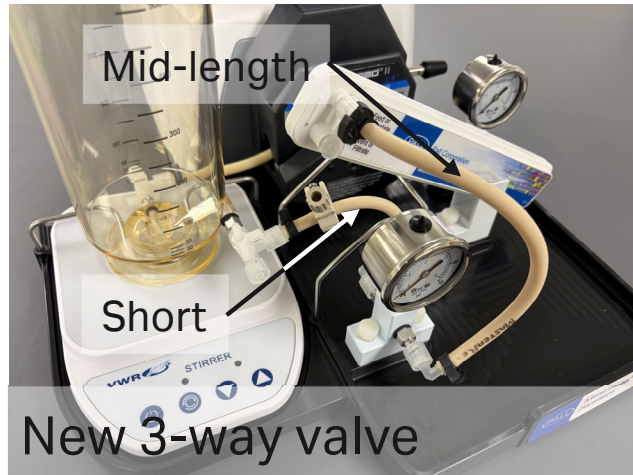
Old connections

Mid-length tube is connected to the rear fitting of the pressure gauge 2 mount and to the retentate port of capsule filter.

Short tube is connected the front fitting of the pressure gauge 2 mount and to port 1 of the 3-way valve on the reservoir inlet.



New connections (to keep the sample port accessible)

Mid-length tube is connected to the front fitting of the pressure gauge 2 mount and to the retentate port of capsule filter.

Short tube is connected the rear fitting of the pressure gauge 2 mount and to port 2 of the 3-way valve on the reservoir inlet

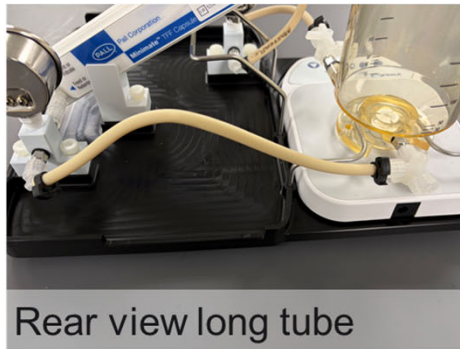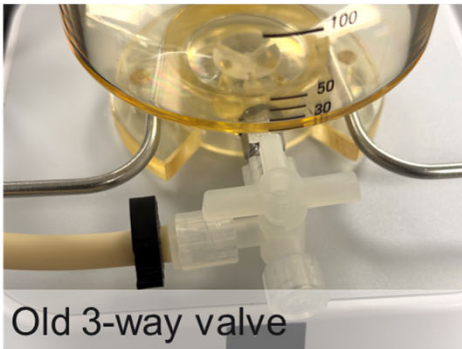Traditional connection of 3-way valve at reservoir rear port

The long tube is connected to port 2 of the 3-way valve at the reservoir outlet via a straight luer lock hose barb and to the rear fitting on the pressure gauge 1 mount via an elbow.


Recommendation for 3-way valve at reservoir rear port – Option 1

The 3-way valve is mounted upside down so port 2 is facing towards the pressure gauge.

The long tube (length 8½") is prepared as customary with male luer lock-1/8" hose barb elbow on one end and a straight male luer lock-1/8" hose barb on the other.

The long tube is connected in the traditional way to port 2 of the 3-way valve at the reservoir outlet via a straight luer lock hose barb and to the rear fitting on the pressure gauge 1 mount via an elbow.

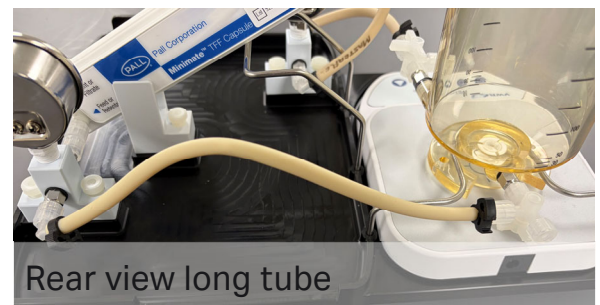This configuration keeps the feed line in the usual position but makes the 3-way valve operation more difficult.

Recommendation for 3-way valve at reservoir rear port – Option 2

The 3-way valve is mounted in normal position with port 2 facing away from the pressure gauge.

The long tube (length 9¼") is prepared with two male luer lock-1/8" hose barb elbows.

The long tube is connected via elbows to both port 1 of the 3-way valve at the reservoir outlet and the rear fitting on the pressure gauge 1 mount.

This configuration causes the feed line to be in a slightly diagonal position. The line will have a slight bend in it when inserted into the pump head (no effect on function) but will allow normal operation of the 3-way valve.

This method requires additional male luer lock-1/8" hose barb elbow that presently is not present in fittings kit or accompanying the system.

Long tube with 2 hose barb elbows

Long tube mounted to rear port

Tubing in pump head – note offset